pH-Fix 0-14



Test strips for the determination of the pH-Value in the range of pH 0-14

Contents:

100 Test strips

1 Colour scale pH 0-14

Instructions for use:

- 1. Dip test strip for 1 sec. with all test fields in the sample solution.
- 2. Shake off excess sample solution.
- 3. Compare with colour scale and read off the corresponding pH-value.

Note:

If sample solution is weakly buffered leave test strip in solution until colour no longer changes.

Storage:

Avoid exposing the test strips to sunlight and moisture. Store the container below + 30 $^{\circ}\text{C}$ in a dry place.

Disposal:

Please dispose used test strips in domestic waste.









MACHEREY-NAGEL GmbH & Co. KG · Neumann-Neander-Str. 6–8 · 52355 Düren · Germany
Tel.: +49 24 21 969-0 · Fax: +49 24 21 969-199 · info@mn-net.com · www.mn-net.com