

## Brix meter Measure 0-55%



### Wide Application :

Refractometer can measure the sugar concentration in fruit, beverage and vegetables, can be used in beverages quality management, food industrial sector, agricultural production scientific research.

It can also be used in the manufacture of beer and wine.

### Specifications :

Temperature compensation:10-40°C

Ambient temperature:10-40°C

Full of electricity:About 20,000 times

Resolution:0.1% Brix

Accuracy:  $\pm 0.2\%$  Brix/ $\pm 1^\circ\text{C}$

Range:0-55%

Battery:2\*AA Battery (cannot be recharged)

Size and weight: 121(L)\*58.5(W)\*30.0(H),119g

# Measurement steps



**01** Power on



**02** Drop the sample, at least 2 / 3 of the sample tank



**03** Press the "meas" key, results in 5 seconds



**04** Clean and dry the sample tank for the next measurement

